PANSERIN™ 401

PANSERIN™ 401 is a complete medium ready for use for the serumfree cultivation of a multitude of cells. Among others the following cells have been cultivated successfully:

- Hybridoma
- Lymphocytes
- Macrophages
- Fibroblasts
- Melanocytes
- Carcinome cells
- HEK-cells
- HeLa-cells
- CHO-cells

Composition:

Based on Iscove's MEM, trace elements, albumin, cholesterol, soya lipids and vitamins were added to the medium. It does not contain any growth or attachment factors or any insulin.

Application:

For many cells (SP2, HEK, L929, CHO) a slow adaption to PANSERIN™ 401 is not necessary.

They are adapted to the serumfree culture according to the **Direct Adaption Method**.

- Transfer vital cells (> 90 % Vitality) of the logarithmic growth phase from the culture containing serum directly into PANSERIN™ 401. A higher cell seed (5×10^4 - 1×10^5) facilitates the adaption to the serumfree culture.
- After 36 - 48 hours in the culture, replace medium by fresh PANSERIN™ 401.
- As soon as the cells have confluent grown at 90 % to 100 %, passage the cells with the usual trypsinating technique. You have to be aware that cells in the serumfree culture show an extremeley sensitive reaction on trypsin. Normal trypsin concentrations (0,25 %) can be used, however the incubation period should be as short as possible at 4 °C. The trypsin can be removed by washing the cells and centrifugation. Soybean-trypsin-inhibitor has to be used with caution as it is toxic for some cells.
- After several passages into PANSERIN™ 401 the adaption is completed.

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PAN™
BIOTECH GmbH

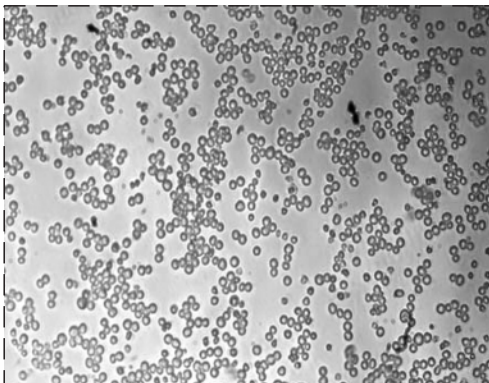
Critical cells can be adapted to the serumfree culture according to the **Gradual Adaption Method**.

- Transfer vital cells (> 90 % Vitality) of the logarithmic growth phase from the culture containing serum (10 % - 20 % FCS) directly into PANSERIN™ 401. Reduce serum contents to 5 %. The usual seeding densities can be kept, but with very critical cells increase seeding densities by 1,5 to 2 x.
- After 2 passages with 5 % FCS supplementation, reduce FCS contents to 1 %.
- After 2 - 3 passages with 1% FCS, reduce serum contents to 0,5 %.
- After further 2 - 3 passages with 0,5 % FCS, cultivate cells without addition of FCS in PANSERIN™ 401.

PANSERIN™ 401 does not contain any attachment factors. With some cell types a pretreatment of the incubation dishes with gelatin, collagen, poly-D-lysine or fibronectin can considerably facilitate the culture under serumfree conditions or even enable it. Please note this above all for low seeding densities.

Additional growth factors are necessary for some cells and have to be added depending on the demand. Please note another important point:

With every adaption to serumfree media, changes of the cells should be taken into consideration. These changes can concern the morphology, the karyotype, the surface marker etc. Thus cells in serumfree medium don't always have to be identical with those from the culture containing serum in which they originate (selection).

SP2/0-Ag-14 in PANSERIN™ 401

Cells were cultivated in PANSERIN™ 401 without adaption

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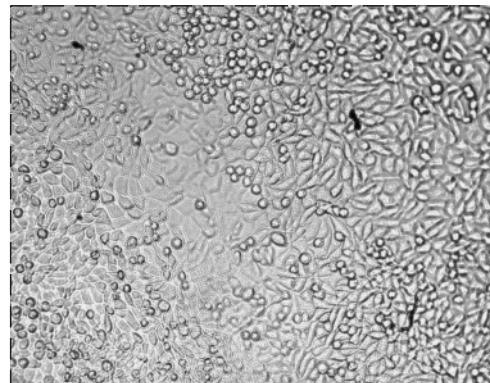
L 929 in PANSERIN™ 401

Cells were cultivated in PANSERIN™ 401 without adaption phase.

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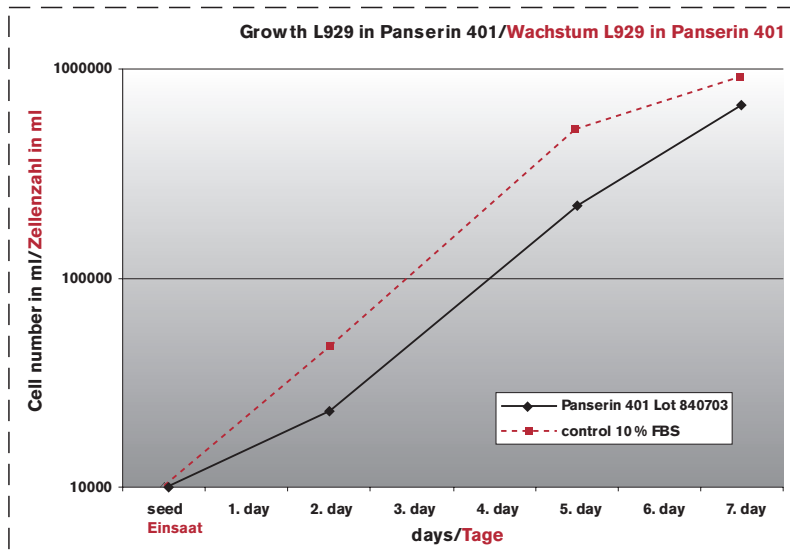
PAN™
BIOTECH GmbH


Typical growth curve L929 in PANSERIN™ 401:

L929 were cultivated in PANSERIN™ 401 without adaption phase. As a comparison L929 in DMEM with 4,5 g/l glucose with 10 % FBS

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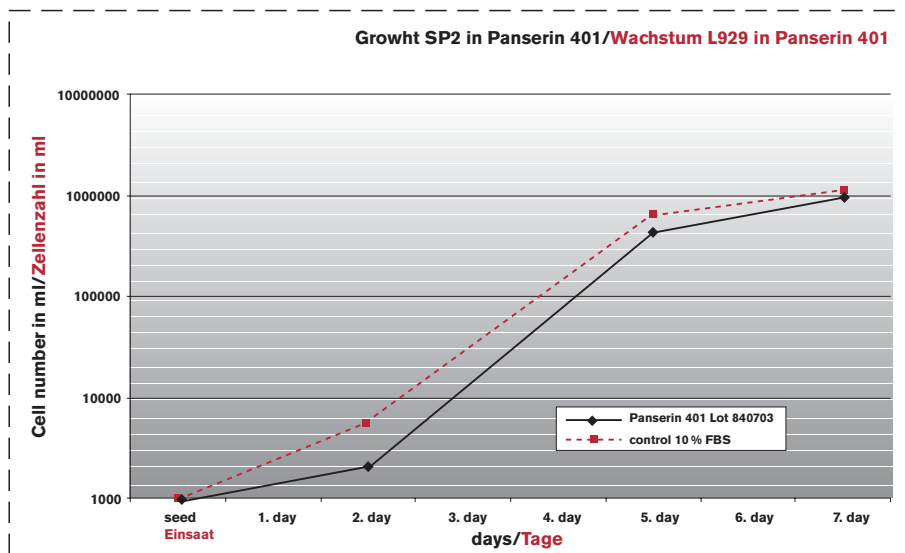Typical growth curve SP2/0-Ag-14 in PANSERIN™ 401:

SP2/0-Ag-14 cells were cultivated in PANSERIN™ 401 without adaption phase. As a comparison cell growth in RPMI 1640 with 10 % FBS

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